



INSTASLIDE SLIDE SHARING

---



## **Microvisioneer InstaSlide - Scan and Share**

A telemicroscopy solution

<https://www.microvisioneer.com/instaslide>



# The Microvisioneer InstaSlide Software


Share digital slides within seconds with anyone, anywhere

Microvisioneer InstaSlide is a software to share folders of virtual slides instantly with remote parties. InstaSlide makes virtual slides available as websites immediately without uploading them to a cloud server.

It is a simple slide sharing solution which works especially well in areas with slow internet connections or in closed local networks.

The slides can be viewed from any distant location, with any device with a browser. Together with the Microvisioneer mvSlide scanning software, InstaSlide represents a complete telemicroscopy solution.

## Slide sharing with InstaSlide

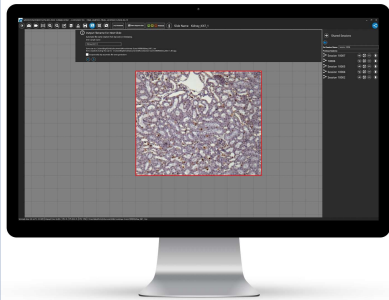


1

**Install and run InstaSlide on your computer.** Start the application. InstaSlide needs to run on the computer that contains the digital slides to be shared. InstaSlide can be fully integrated into Microvisioneer's manual scanning software as scans are directly saved into predefined sessions folders.

2

**Select and share scans.** For each scanning session, a worldwide accessible URL is created within seconds. Select the contacts supposed to view the slides in your personalized address book. The selected contacts automatically receive an e-mail notification within seconds containing the URL.

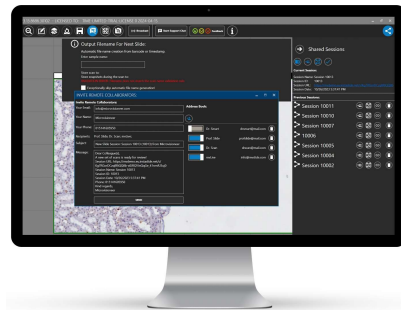


3

**View the shared zoomable whole slide images.** Everyone receiving the URL can view the virtual slides now. Nothing else than a web browser is needed. The scans can be viewed as a zoomable website on any device such as computers, tablets or smartphones. The capacity of the InstaSlide slide server is unlimited. You can share hundreds of slides within seconds.



# The Digital Workflow with InstaSlide

The easy transition to fully digital workflows

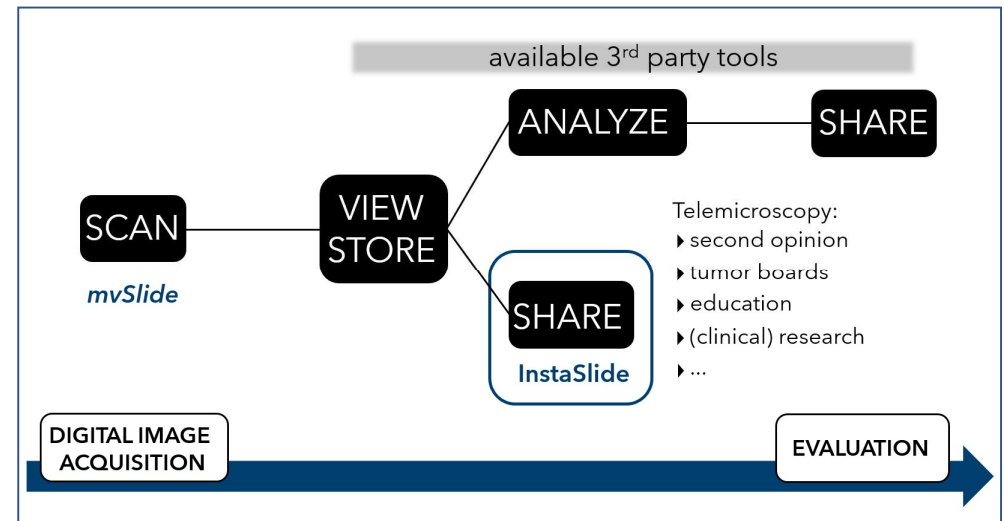

## Benefits of InstaSlide



### Read slides anywhere

There is no need to visit remote sites. Save your precious time and simply ask a colleague to scan slides and share them instantly via InstaSlide. Slides are accessible with any device at any location.



### Share slides with anyone

Share difficult slides with your colleagues for a second opinion. They just need a web browser to have a look at the scans.



### Save time

Common upload of scans to cloud platforms takes a lot of time. InstaSlide shares even terabytes of virtual slides within seconds. Sharing and reading slides via InstaSlide is extremely fast, even with slow internet connections.



### Maintain image quality

The InstaSlide web viewer does not use the traditional screen sharing technology often used by telemicroscopy solutions. As a result, the image quality is superior to traditional techniques.

# Specifications


Key technical features of Microvisioneer InstaSlide

InstaSlide Versions	Microvisioneer InstaSlide makes folders of virtual slides available as websites without uploading them to a cloud server. There are two versions of InstaSlide available.
Standard Version	Makes scans available via a web URL
LAN Edition	Makes scans available in a local network only
Supported file types	InstaSlide supports sharing images in *.svs and *.tiff file format
Scan accessibility	Scans are accessible while InstaSlide is running. Disable automatic sleep mode of computer; Add InstaSlide to autostart configuration
Supported remote devices	Any device with a browser
InstaSlide URL validity	Infinite, also after computer restart
InstaSlide URL encryption	Unique, encrypted URL creation
Possibility of integration into Microvisioneer's slide scanning software	InstaSlide can be directly integrated into mvSlide for easy and immediate slide sharing via a personalized address book and automatically created e-mail notifications.

Microvisioneer InstaSlide is for research use only.


# Microvisioneer Product Range

Find your solution

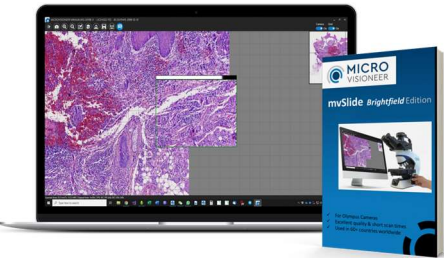

## Microvisioneer mvSlide Brightfield Edition

Manual scanning for brightfield applications, oil immersion, polarized light and phase contrast. Suitable for histology, immunohistochemistry, cytology, material science and many more.

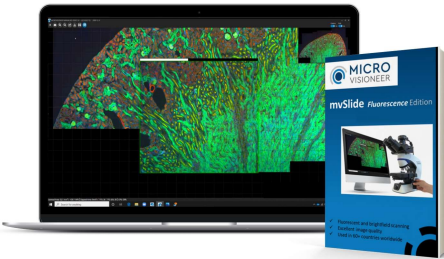

## Microvisioneer mvSlide Fluorescence Edition

Manual scanning for *brightfield and fluorescence* applications. Both single channel and multi channel fluorescence imaging is possible.



## Microvisioneer InstaSlide Slide Sharing Software

Immediate sharing of virtual slides via the internet with anyone, anywhere. A Microvisioneer telemicroscopy solution.




## Microvisioneer mvLive Broadcasting Software

Live sharing of the camera image with remote colleagues. Best image quality without compression artefacts.

**Microvisioneer GmbH**  
Germany  
VAT ID: DE366777766

---

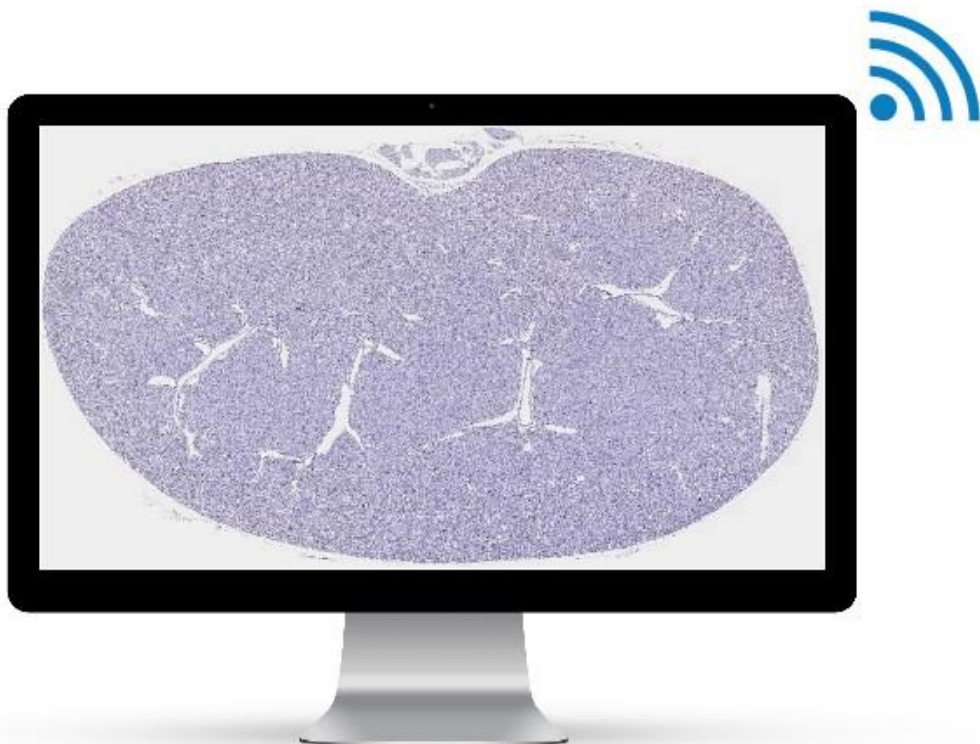

Contact us to discuss your opportunities or to book an online demo.

Email [info@microvisioneer.com](mailto:info@microvisioneer.com)  
Phone +49(0) 151 41605950

<https://www.microvisioneer.com>

